



碧云天网站



微信公众号

碧云天生物技术/Beyotime Biotechnology

订货热线: 400-168-3301或800-8283301

订货e-mail: order@beyotime.com

技术咨询: info@beyotime.com

网址: http://www.beyotime.com

## Cy3标记山羊抗兔IgG (H+L)

产品编号	产品名称	包装
A0516	Cy3标记山羊抗兔IgG (H+L)	100 $\mu$ l

### 产品简介:

- 本Cy3标记山羊抗兔IgG (H+L) (Cy3-labeled Goat Anti-Rabbit IgG (H+L))为进口分装, 用于免疫荧光染色。
- Cy3是近年新开发的一种红色荧光探针。它比绝大部分其它的红色荧光探针更加明亮, 更加不容易淬灭, 而且背景更低。Cy3的吸收(激发)和发射峰参见下表。

Fluorophore	Absorption Peak (nm)	Emission Peak (nm)
Cy3	550	570

- 本抗体为用纯化的兔IgG免疫山羊, 然后用亲和纯化柱对获得的抗血清进行纯化, 并经过人IgG吸附纯化的优质二抗。对人IgG几乎没有结合能力。
- 本Cy3标记山羊抗兔IgG (H+L)用于免疫荧光染色时的推荐稀释比例为1:500。实际实验操作过程中需根据抗原和抗体的具体情况适当调节荧光标记二抗的稀释比例, 推荐的调节范围为1:200-1000。
- 本抗体如果用于常规的免疫染色, 以每次检测需1毫升1:500稀释的荧光标记二抗计, 至少可以检测50次。如果适当重复使用已经使用过的荧光标记二抗, 至少可以多检测150-250次。

### 包装清单:

产品编号	产品名称	包装
A0516	Cy3标记山羊抗兔IgG (H+L)	100 $\mu$ l
—	说明书	1份

### 保存条件:

-20 $^{\circ}$ C避光保存, 一年有效。

### 注意事项:

- 本产品仅限于专业人员的科学研究用, 不得用于临床诊断或治疗, 不得用于食品或药品, 不得存放于普通住宅内。
- 为了您的安全和健康, 请穿实验服并戴一次性手套操作。

### 使用说明:

1. 免疫荧光染色请参考相关实验步骤进行。起始稀释浓度按照产品简介中推荐的稀释比例进行稀释。
2. 如果希望重复使用稀释的荧光标记二抗, 稀释的荧光标记二抗4 $^{\circ}$ C保存。

### 使用本产品的文献:

1. Liao H, Huang W, Schachner M, Guan Y, Guo J, Yan J, Qin J, Bai X, Zhang L. Beta 1 integrin-mediated effects of tenascin-R domains EGFL and FN6-8 on neural stem/progenitor cell proliferation and differentiation in vitro. *J Biol Chem.* 2008 Oct 10;283(41):27927-36.
2. Zhu J, Zhou Z, Liu Y, Zheng J. Fractalkine and CX3CR1 are involved in the migration of intravenously grafted human bone marrow stromal cells toward ischemic brain lesion in rats. *Brain Res.* 2009 Sep 1;1287:173-83.
3. Li W, Cun W, Liu L, Hong M, Wang L, Wang L, Dong C, Li Q. The transactivating effect of HSV-1 ICP0 is enhanced by its interaction with the PCAF component of histone acetyltransferase. *Arch Virol.* 2009;154(11):1755-64.
4. Wong I, Liao H, Bai X, Zaknic A, Zhong J, Guan Y, Li HY, Wang YJ, Zhou XF. ProBDNF inhibits infiltration of ED1+ macrophages after spinal cord injury. *Brain Behav Immun.* 2010;24(4):585-97.
5. Chen J, Song M, Yu S, Gao P, Yu Y, Wang H, Huang L. Advanced glycation end products alter functions and promote apoptosis in endothelial progenitor cells through receptor for advanced glycation end products mediate overexpression of cell oxidant stress. *Mol Cell Biochem.* 2010;335(1-2):137-46.
6. Zhang Y, Zhou L, Bao YL, Wu Y, Yu CL, Huang YX, Sun Y, Zheng LH, Li YX. Butyrate induces cell apoptosis through activation of JNK MAP kinase pathway in human colon cancer RKO cells. *Chem Biol Interact.* 2010;185(3):174-81.
7. Hu B, Yang Y, Liu J, Ma Z, Huang H, Liu S, Yu Y, Hao Y, Wang B, Lu M, Yang D. Establishment of a functional cell line expressing both subunits of H1a and H2c of human hepatocyte surface molecule ASGPR. *J Huazhong Univ Sci Technol Med Sci.* 2010 Oct;30(5):556-61.
8. Zheng Y, Zhou M, Ye A, Li Q, Bai Y, Zhang Q. The conformation change of Bcl-2 is involved in arsenic trioxide-induced apoptosis and inhibition of proliferation in SGC7901 human gastric cancer cells. *World J Surg Oncol.* 2010 Apr 20;8:31.
9. Li W, Wang G, Zhang H, Xin G, Zhang D, Zeng J, Chen X, Xu Y, Cui Y, Li K. Effects of NS1 variants of H5N1 influenza virus on interferon induction, TNFalpha response and p53 activity. *Cell Mol Immunol.* 2010 May;7(3):235-42.
10. Zhang Y, Bao YL, Yang MT, Wu Y, Yu CL, Huang YX, Sun Y, Zheng LH, Li YX. Activin A induces SLC5A8 expression through the Smad3 signaling pathway in human colon cancer RKO cells. *Int J Biochem Cell Biol.* 2010 Dec;42(12):1964-72.
11. Zhang Y, Zhou L, Bao YL, Wu Y, Yu CL, Huang YX, Sun Y, Zheng LH, Li YX. Butyrate induces cell apoptosis through activation of JNK MAP kinase pathway in human colon cancer RKO cells. *Chem Biol Interact.* 2010 May 14;185(3):174-81.
12. Meng G, Liu Y, Lou C, Yang H. Emodin suppresses lipopoly saccharide-induced pro-inflammatory responses and NF- $\kappa$ B activation by disrupting lipid rafts in CD14-negative endothelial cells. *Br J Pharmacol.*

13. 王福科, 刘流, 赵德萍, 何晓光, 代晓明, 李逸松, 李玉晓. 体外血管内皮细胞和脂肪干细胞联合培养对脂肪干细胞成骨分化的影响. 中国修复重建外科. 2010;Vol.14(32):6001-5.
14. Zheng Y, Zhou M, Ye A, Li Q, Bai Y, Zhang Q. The conformation change of Bcl-2 is involved in arsenic trioxide-induced apoptosis and inhibition of proliferation in SGC7901 human gastric cancer cells. *World J Surg Oncol.* 2010 Apr 20;8:31.
15. Wang F, Wang F, Zou Z, Liu D, Wang J, Su Y. Active deformation of apoptotic intestinal epithelial cells with adhesion-restricted polarity contributes to apoptotic clearance. *Lab Invest.* 2011 Mar;91(3):462-71.
16. 景黎君, 贾永林, 鲁晶晶, 韩瑞, 王舒阳, 李尽义. MicroRNA-9-1慢病毒载体的构建及其对小鼠骨髓间质干细胞诱导分化为神经细胞的影响. 中国病理生理杂志. 2011;Vol.27(2):326-331.
17. Yang L, Ping YF, Yu X, Qian F, Guo ZJ, Qian C, Cui YH, Bian XW. Gastric cancer stem-like cells possess higher capability of invasion and metastasis in association with a mesenchymal transition phenotype. *Cancer Lett.* 2011 Nov 1;310(1):46-52.
18. Yu PJ, Ci W, Wang GF, Zhang JY, Wu SY, Xu W, Jin H, Zhu ZG, Zhang JJ, Pang JX, Wu SG. Praeruptorin A inhibits lipopolysaccharide-induced inflammatory response in murine macrophages through inhibition of NF- $\kappa$ B pathway activation. *Phytother Res.* 2011 Apr;25(4):550-6.
19. Yu PJ, Jin H, Zhang JY, Wang GF, Li JR, Zhu ZG, Tian YX, Wu SY, Xu W, Zhang JJ, Wu SG. Pyranocoumarins isolated from *Peucedanum praeruptorum* Dunn suppress lipopolysaccharide-induced inflammatory response in murine macrophages through inhibition of NF- $\kappa$ B and STAT3 activation. *Inflammation.* 2012 Jun;35(3):967-77.
20. Jin H, Zhu ZG, Yu PJ, Wang GF, Zhang JY, Li JR, Ai RT, Li ZH, Tian YX, Zhang WX, Wu SG. Myrislignan Attenuates Lipopoly saccharide-induced Inflammation Reaction in Murine Macrophage Cells Through Inhibition of NF- $\kappa$ B Signalling Pathway Activation. *Phytother Res.* 2012 Sep;26(9):1320-6.
21. Huo SJ, Li YC, Xie J, Li Y, Raisman G, Zeng YX, He JR, Weng CH, Yin ZQ. Transplanted Olfactory Ensheathing Cells Reduce Retinal Degeneration in Royal College of Surgeons Rats. *Curr Eye Res.* 2012 Aug;37(8):749-58.
22. Huang S, Lu G, Wu Y, Jirigala E, Xu Y, Ma K, Fu X. Mesenchymal stem cells delivered in a microsphere-based engineered skin contribute to cutaneous wound healing and sweat gland repair. *J Dermatol Sci.* 2012 Apr;66(1):29-36.
23. Zhu L, Yuan H, Guo C, Lu Y, Deng S, Yang Y, Wei Q, Wen L, He Z. Zearalenone induces apoptosis and necrosis in porcine granulosa cells via a caspase-3- and caspase-9-dependent mitochondrial signaling pathway. *J Cell Physiol.* 2012 May;227(5):1814-20.
24. Liu Y, Yu J, Du M, Wang W, Zhang W, Wang Z, Jiang X. Accelerating microfluidic immunoassays on filter membranes by applying vacuum. *Biomed Microdevices.* 2012 Feb;14(1):17-23.
25. Guo HF, Li ZS, Dong SW, Chen WJ, Deng L, Wang YF, Ying DJ. Piezoelectric PU/PVDF electrospun scaffolds for wound healing applications. *Colloids Surf B Biointerfaces.* 2012 Aug 1;96:29-36.
26. Xue H, Zhou L, Yuan P, Wang Z, Ni J, Yao T, Wang J, Huang Y, Yu C, Lu L. Counteraction between angiotensin II and angiotensin-(1-7) via activating angiotensin type I and Mas receptor on rat renal mesangial cells. *Regul Pept.* 2012 Aug 20;177(1-3):12-20.
27. Li Y, Li C, Chen Z, He J, Tao Z, Yin ZQ. A microRNA, mir133b, suppresses melanopsin expression mediated by failure dopaminergic amacrine cells in RCS rats. *Cell Signal.* 2012 Mar;24(3):685-98.
28. Zhang HJ, Yao DF, Yao M, Huang H, Wang L, Yan MJ, Yan XD, Gu X, Wu W, Lu SL. Annexin A2 silencing inhibits invasion, migration, and tumorigenic potential of hepatoma cells. *World J Gastroenterol.* 2013 Jun 28;19(24):3792-801.
29. Zhang B, Cui S, Bai X, Zhuo L, Sun X, Hong Q, Fu B, Wang J, Chen X, Cai G. SIRT3 overexpression antagonizes high glucose accelerated cellular senescence in human diploid fibroblasts via the SIRT3-FOXO1 signaling pathway. *Age (Dordr).* 2013 Dec;35(6):2237-53.
30. Li X, Li Z, Li N, Qi J, Fan K, Yin P, Zhao C, Liu Y, Yao W, Cai X, Wang L, Zha X. MAGI2 enhances the sensitivity of BEL-7404 human hepatocellular carcinoma cells to staurosporine-induced apoptosis by increasing PTEN stability. *Int J Mol Med.* 2013 Aug;32(2):439-47.
31. Li Q, Liu QH, Huang J. F0ATP synthase b-chain of *Litopenaeus vannamei* involved in White Spot Syndrome Virus infection. 2013 Aug;47(1):42-8.
32. Liu H, Xu H, Yu T, Yao J, Zhao C, Yin ZQ. Expression of Perineuronal Nets, Parvalbumin and Protein Tyrosine Phosphatase  $\sigma$  in the Rat Visual Cortex During Development and After BFD. *Curr Eye Res.* 2013 Oct;38(10):1083-94.
33. Zhou S, Tang L, Wang H, Dai J, Zhang J, Shen L, Ng SW, Berkowitz RS. Overexpression of c-Abl predicts unfavorable outcome in epithelial ovarian cancer. *Gynecol Oncol.* 2013 Oct;131(1):69-76.
34. Li Q, Zhang Z, Li Z, Zhou M, Liu B, Pan L, Ma Z, Zheng Y. ADAM17 Is Critical for Multipolar Exit and Radial Migration of Neuronal Intermediate Progenitor Cells in Mice Cerebral Cortex. *PLoS One.* 2013 Jun 3;8(6):e65703.
35. Yang N, Hui L, Wang Y, Yang H, Jiang X. Overexpression of SOX2 promotes migration, invasion, and epithelial-mesenchymal transition through the Wnt/ $\beta$ -catenin pathway in laryngeal cancer Hep-2 cells. *Tumour Biol.* 2014 Aug;35(8):7965-73.
36. Peng H, Zhang W, Xiao T, Zhang Y. Nlrp4g is an oocyte-specific gene but is not required for oocyte maturation in the mouse. *Reprod Fertil Dev.* 2014 Jun;26(5):758-68.
37. Li H, Zhang X, Zeng S, Chen L, Li X, Lin C, Zhang M, Shu S, Xie S, He Y, Yang L, Tang S, Fu X. The cellular localization of Na<sup>+</sup>/H<sup>+</sup> exchanger 1, cystic fibrosis transmembrane conductance regulator, potassium channel, epithelial sodium channel  $\gamma$  and vacuolar-type H<sup>+</sup>-ATPase in human eccrine sweat glands. *Acta Histochem.* 2014 Oct;116(8):1237-43.
38. Zhao F, Li L, Guan L, Yang H, Wu C, Liu Y. Roles for GP IIb/IIIa and  $\alpha$ v $\beta$ 3 integrins in MDA-MB-231 cell invasion and shear flow-induced cancer cell mechanotransduction. *Cancer Lett.* 2014 Mar 1;344(1):62-73.
39. Zhang JR, Yao FH, Zhang JG, Ji ZY, Li KL, Zhan J, Tong YN, Lin LR, He YN. Ischemia-reperfusion induces renal tubule pyroptosis via the CHOP-caspase-11 pathway. *Am J Physiol Renal Physiol.* 2014 Jan 1;306(1):F75-84.
40. Li H, Chen L, Zeng S, Li X, Zhang X, Lin C, Zhang M, Xie S, He Y, Shu S, Yang L, Tang S, Fu X. Matrigel basement membrane matrix induces eccrine sweat gland cells to reconstitute sweat gland-like structures in nude mice. *Exp Cell Res.* 2015 Mar 1;332(1):67-77.
41. Zhang L, Li H2, Zeng S, Chen L, Fang Z, Huang Q. Long-term tracing of the BrdU label-retaining cells in adult rat brain. *Neurosci Lett.* 2015 Mar 30;591:30-4.
42. Liang GJ, Zhang XF, Wang JJ, Sun YC, Sun XF, Cheng SF, Li L, De Felici M, Shen W. Activin A Accelerates the Progression of Fetal Oocytes Throughout Meiosis and Early Oogenesis in the Mouse. *Stem Cells Dev.* 2015 Oct 15;24(20):2455-65.
43. Si J, Dai J, Zhang J, Liu S, Gu J, Shi J, Shen SG, Guo L. Comparative investigation of human amniotic epithelial cells and mesenchymal stem cells for application in bone tissue engineering. *Stem Cells Int.* 2015;2015:565732.
44. Wang QZ, Gao HQ, Liang Y, Zhang J, Wang J, Qiu J. Cofilin1 is involved in hypertension-induced renal damage via the regulation of NF- $\kappa$ B in renal tubular epithelial cells. *J Transl Med.* 2015 Oct 8;13(1):323.
45. Jin Y, Lin L, Lin Y, Zheng C. Exogenous carcinoembryonic antigen-related cell adhesion molecule 1 suppresses 2,4,6-trinitrobenzene sulfonic acid-induced ulcerative colitis in mice. *J Surg Res.* 2015 May 1;195(1):113-20.
46. Zhan Y, Xiang F, Wu R, Xu J, Ni Z, Jiang J, Kang X. MiRNA-149 modulates chemosensitivity of ovarian cancer A2780 cells to paclitaxel by targeting MyD88. *J Ovarian Res.* 2015 Jul 30;8:48.
47. Cao Q, Jiang Y, Shi J, Xu C, Liu X, Yang T, Fu P, Niu T. Artemisinin inhibits the proliferation, migration, and inflammatory reaction induced by tumor necrosis factor- $\alpha$  in vascular smooth muscle cells through nuclear factor kappa B pathway. *J Surg Res.* 2015 Apr;194(2):667-78.
48. Dai C, Tang S, Velkov T, Xiao X. Colistin-Induced Apoptosis of Neuroblastoma-2a Cells Involves the Generation of Reactive Oxygen Species, Mitochondrial Dysfunction, and Autophagy. *Mol Neurobiol.* 2016 Sep;53(7):4685-700.
49. Xia X, Feng H, Li C, Qin C, Song Y, Zhang Y, Lan X. <sup>99m</sup>Tc-labeled estradiol as an estrogen receptor probe: Preparation and preclinical evaluation. *Nucl Med Biol.* 2016 Jan;43(1):89-96.
50. Li H, Chen L, Zhang M, Zhang B. Foxa1 gene and protein in developing rat eccrine sweat glands. *J Mol Histol.* 2016 Oct 27. [Epub ahead of print]
51. Li H, Li X, Zhang B, Zhang M, Chen W, Tang S, Fu X. Changes in keratins and alpha-smooth muscle actin during three-dimensional reconstitution of eccrine sweat glands. *Cell Tissue Res.* 2016

Jul;365(1):113-22.

52. Hou C, Wu Q, Ouyang C, Huang T. Effects of an intravitreal injection of interleukin-35-expressing plasmid on pro-inflammatory and anti-inflammatory cytokines. *Int J Mol Med*. 2016 Sep;38(3):713-20.
53. Dong Z, Song Q, Zhang Y, Chen S, Zhang X, Zhao P, Xia Q. Structure, evolution, and expression of antimicrobial silk proteins, seroins in Lepidoptera. *Insect Biochem Mol Biol*. 2016 Aug;75:24-31.
54. Li H, Zhang M, Chen L, Li X, Zhang B. Human eccrine sweat gland cells reconstitute polarized spheroids when subcutaneously implanted with Matrigel in nude mice. *J Mol Histol*. 2016 Oct;47(5):485-90.
55. Jin F, Qiao C, Luan N, Li H. Lentivirus-mediated PHLDA2 overexpression inhibits trophoblast proliferation, migration and invasion, and induces apoptosis. *Int J Mol Med*. 2016 Apr;37(4):949-57.
56. Zhou Y, Zhang S, Dai C, Tang S, Yang X, Li D, Zhao K, Xiao X. Quinocetone triggered ER stress-induced autophagy via ATF6/DAPK1-modulated mAtg9a trafficking. *Cell Biol Toxicol*. 2016 Apr;32(2):141-52.
57. Zhou Y, He Z, Gao Y, Zheng R, Zhang X, Zhao L, Tan M. Induced Pluripotent Stem Cells Inhibit Bleomycin-Induced Pulmonary Fibrosis in Mice through Suppressing TGF- $\beta$ 1/Smad-Mediated Epithelial to Mesenchymal Transition. *Front Pharmacol*. 2016 Nov 15;7:430.
58. Gao Z, Han Y, Hu Y, Wu X, Wang Y, Zhang X, Fu J, Zou X, Zhang J, Chen X, Jose PA, Lu X, Zeng C. Targeting HO-1 by Epigallocatechin-3-Gallate Reduces Contrast-Induced Renal Injury via Anti-Oxidative Stress and Anti-Inflammation Pathways. *PLoS One*. 2016 Feb 11;11(2):e0149032.
59. Xia X, Feng H, Li C, Qin C, Song Y, Zhang Y, Lan X. <sup>99m</sup>Tc-labeled estradiol as an estrogen receptor probe: Preparation and preclinical evaluation. *Nucl Med Biol*. 2016 Jan;43(1):89-96.
60. Zhang H, Ju B, Zhang X, Zhu Y, Nie Y, Xu Y, Lei Q. Magnolol Attenuates Concanavalin A-induced Hepatic Fibrosis, Inhibits CD4+ T Helper 17 (Th17) Cell Differentiation and Suppresses Hepatic Stellate Cell Activation: Blockade of Smad3/Smad4 Signaling. *Basic Clin Pharmacol Toxicol*. 2016 Dec 29.
61. Zhang H, Nan W, Wang S, Zhang T, Si H, Yang F, Li G. Epidermal Growth Factor Promotes Proliferation and Migration of Follicular Outer Root Sheath Cells via Wnt/ $\beta$ -Catenin Signaling. *Cell Physiol Biochem*. 2016;39(1):360-70.
62. Dai C, Ciccotosto GD, Cappai R, Tang S, Li D, Xie S, Xiao X, Velkov T. Curcumin Attenuates Colistin-Induced Neurotoxicity in N2a Cells via Anti-inflammatory Activity, Suppression of Oxidative Stress, and Apoptosis. *Mol Neurobiol*. 2016 Dec 12. [Epub ahead of print]

Version 2016.10.10