
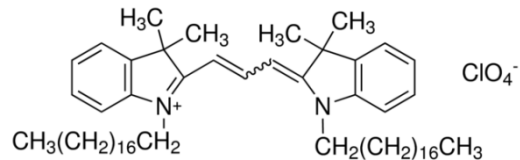


DiI (细胞膜红色荧光探针)

产品编号	产品名称	包装
C1036	DiI (细胞膜红色荧光探针)	10mg

产品简介:

- DiI即DiIC₁₈(3), 全称为1,1'-dioctadecyl-3,3,3',3'-tetramethylindocarbocyanine perchlorate, 是最常用的细胞膜荧光探针之一, 呈现橙红色荧光。DiI是一种亲脂性膜染料, 进入细胞膜后可以侧向扩散逐渐使整个细胞的细胞膜被染色。
- DiI在进入细胞膜之前荧光非常弱, 仅当进入到细胞膜后才可以被激发出很强的荧光。DiI被激发后可以发出橙红色的荧光, DiI和磷脂双层膜结合后的激发光谱和发射光谱参考下图。其中, 最大激发波长为549nm, 最大发射波长为565nm。

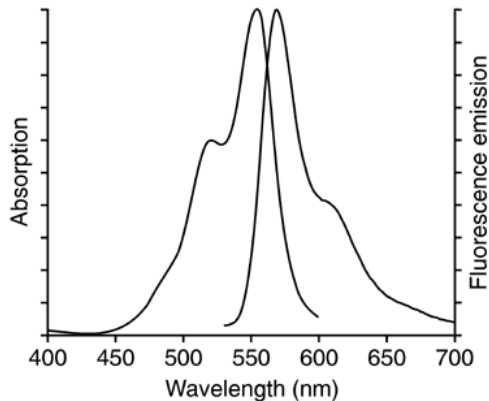

- DiI的分子式为C₅₉H₉₇ClN₂O₄, 分子量为933.88, CAS number为41085-99-8。
- DiI可以溶解于无水乙醇、DMSO和DMF, 其中在DMSO中的溶解度大于10mg/ml。发现较难溶解时可以适当加热, 并用超声处理以促进溶解。
- DiI 被广泛用于正向或逆向的, 活的或固定的神经等细胞或组织的示踪剂或长期示踪剂(long-term tracer)。DiI通常不会影响细胞的生存力(viability)。被DiI标记的神经细胞在体外培养的条件下可以存活长达4周, 在体内可以长达一年。DiI在经过固定的神经元细胞膜上的迁移速率为0.2-0.6mm/day, 在活的神经元细胞膜上的迁移速率为6mm/day。
- DiI除了最简单的细胞膜荧光标记外, 还可以用于检测细胞的融合和粘附, 检测发育或移植过程中细胞迁移, 通过FRAP(Fluorescence Recovery After Photobleaching)检测脂在细胞膜上的扩散, 检测细胞毒性和标记脂蛋白等。
- 用于细胞膜荧光标记时, DiI的常用浓度为1-25μM, 最常用的浓度为5-10μM。DiI可以直接染色活的细胞或组织, 染色时间通常为5-20分钟。对于固定的细胞或组织, 通常宜使用配制在PBS中的4%多聚甲醛进行固定, 使用其它不适当的固定液会导致荧光背景较高。

包装清单:

产品编号	产品名称	包装
C1036	DiI (细胞膜红色荧光探针)	10mg
—	说明书	1份

保存条件:

4°C避光保存, 一年有效。配制的储存液-20°C避光保存, 半年有效。

注意事项:

- 荧光染料均存在淬灭问题, 请尽量注意避光, 以减缓荧光淬灭。
- 本产品仅限于专业人员的科学研究用, 不得用于临床诊断或治疗, 不得用于食品或药品, 不得存放于普通住宅内。
- 为了您的安全和健康, 请穿实验服并戴一次性手套操作。

使用本产品的文献:

- Zhu C, Huang H, Hua R, Li G, Yang D, Luo J, Zhang C, Shi L, Benovic JL, Zhou N. Molecular and functional characterization of adipokinetic hormone receptor and its peptide ligands in *Bombyx mori*. FEBS Lett. 2009 May 6;583(9):1463-8.
- Danhier F, Lecouturier N, Vroman B, Jérôme C, Marchand-Brynaert J, Feron O, Préat V. Paclitaxel-loaded PEGylated PLGA-based nano particles: in vitro and in vivo evaluation. J Control Release. 2009 Jan 5; 133(1):11-7.
- Chen DQ, Jiang XQ, Zhang C, Ping QE. pH-Sensitive mPEG-Hz-Cholesterol Conjugates as a Liposome Delivery System. Journal of

4. Yu L, Lv JC, Zhou XJ, Zhu L, Hou P, Zhang H. Abnormal expression and dysfunction of novel SGLT2 mutations identified in familial renalglucosuria patients. *Hum Genet.* 2011 Mar;129(3):335-44.
5. Li RJ, Ying X, Zhang Y, Ju RJ, Wang XX, Yao HJ, Men Y, Tian W, Yu Y, Zhang L, Huang RJ, Lu WL. All-trans retinoic acid stealth liposomes prevent the relapse of breast cancer arising from the cancer stem cells. *J Control Release.* 2011 Feb 10;149(3):281-91.
6. Liu M, Li ZH, Xu FJ, Lai LH, Wang QQ, Tang GP, Yang WT. An oligopeptide ligand-mediated therapeutic gene nanocomplex for liver cancer-targeted therapy. *Biomaterials.* 2012 Mar;33(7):2240-50.
7. Cheng C, Zhang XG, Xiang JX, Wang YX, Zhang C, Lu ZT, Li CX. Development of novel self-assembled poly(3-acrylamidophenylboronic acid)/poly(2-lactobionamidoethyl methacrylate) hybrid nanoparticles for improving nasal adsorption of insulin. *Soft Matter.* 2012 8;3:765-73.
8. Cheng C, Zhang XG, Wang YX, Sun L, Li CX. Phenylboronic acid-containing block copolymers: synthesis, self-assembly, and application for intracellular delivery of proteins. *New J. Chem.* 2012 Apr;6:1413-21.
9. Zhao W, Zhang WP, Zhang ZL, He RL, Lin Y, Xie M, Wang HZ, Pang DW. Robust and highly sensitive fluorescence approach for point-of-care virus detection based on immunomagnetic separation. *Anal Chem.* 2012 Mar 6;84(5):2358-65.
10. Li G, Lin DH, Xie XX, Qin LF, Wang JT. Cellular Uptake of PEGylated PLGA Nanoparticles in HeLa Cells. *Frontier of Nanoscience and Technology II.* 2012 Jun;(528):80-3.
11. Li G, Zhou Q, Yu Y, Chen L, Shi Y, Luo J, Benovic J, Lu J, Zhou N. Identification and characterization of distinct C-terminal domains of the human hydroxycarboxylic acid receptor-2 that are essential for receptor export, constitutive activity, desensitization, and internalization. *Mol Pharmacol.* 2012 Dec;82(6):1150-61.
12. Bo Y, Yan L, Gang Z, Tao L, Yinghui T. Effect of calcitonin gene-related peptide on osteoblast differentiation in an osteoblast and endothelial cell co-culture system. *Cell Biol Int.* 2012 Oct 1;36(10):909-15.
13. Chen C, Diao D, Guo L, Shi M, Gao J, Hu M, Yu M, Qian L, Guo N. All-trans-retinoic acid modulates ICAM-1 N-glycan composition by influencing GnT-III levels and inhibits cell adhesion and trans-endothelial migration. *PLoS One.* 2012;7(12):e52975.
14. Guo T, Ke L, Qi B, Wan J, Ge J, Bai L, Cheng B. PTX3 is located at the membrane of late apoptotic macrophages and mediates the phagocytosis of macrophages. *J Clin Immunol.* 2012 Apr;32(2):330-9.
15. Yin T, Wang P, Zheng R, Zheng B, Cheng D, Zhang X, Shuai X. Nanobubbles for enhanced ultrasound imaging of tumors. *Int J Nanomedicine.* 2012;7:895-904.
16. Liu L, Kou P, Zeng Q, Pei G, Li Y, Liang H, Xu G, Chen S. CD4+ T Lymphocytes, especially Th2 cells, contribute to the progress of renal fibrosis. *Am J Nephrol.* 2012;36(4):386-96.
17. Meng Z, Zhang XY, Guo J, Xiang LX, Shao JZ. Scavenger receptor in fish is a lipopolysaccharide recognition molecule involved in negative regulation of NF- κ B activation by competing with TNF receptor-associated factor 2 recruitment into the TNF- α signaling pathway. *J Immunol.* 2012 Oct 15;189(8):4024-39.
18. Guo HF, Li ZS, Dong SW, Chen WJ, Deng L, Wang YF, Ying DJ. Piezoelectric PU/PVDF electrospun scaffolds for wound healing applications. *Colloids Surf B Biointerfaces.* 2012 Aug 1;96:29-36.
19. Jiang H, Chen SS, Yang J, Chen J, He B, Zhu LH, Wang L. CREB-binding protein silencing inhibits thrombin-induced endothelial progenitor cells angiogenesis. *Mol Biol Rep.* 2012 Mar;39(3):2773-9.
20. Bi RY, Kou XX, Meng Z, Wang XD, Ding Y, Gan YH. Involvement of trigeminal ganglionic Na(v) 1.7 in hyperalgesia of inflamed temporomandibular joint is dependent on ERK1/2 phosphorylation of glial cells in rats. *Eur J Pain.* 2013 Aug;17(7):983-94.
21. He B, Jia Z, Du W, Yu C, Fan Y, Dai W, Yuan L, Zhang H, Wang X, Wang J, Zhang X, Zhang Q. The transport pathways of polymer nanoparticles in MDCK epithelial cells. *Biomaterials.* 2013 Jun;34(17):4309-26.
22. Li J, Cheng Y, Zhang X, Zheng L, Han Z, Li P, Xiao Y, Zhang Q, Wang F. The in vivo immunomodulatory and synergistic anti-tumor activity of thymosin α 1-thymopentin fusion peptide and its binding to TLR2. *Cancer Lett.* 2013 Sep 1;337(2):237-47.
23. Zhu H, Cao J, Cui S, Qian Z, Gu Y. Enhanced tumor targeting and antitumor efficacy via hydroxycamptothecin-encapsulated folate-modified N-succinyl-N'-octyl chitosan micelles. *J Pharm Sci.* 2013 Apr;102(4):1318-32.
24. Zhang XY, Xiao YQ, Zhang Y, Ye W. Protective effect of pioglitazone on retinal ischemia/reperfusion injury in rats. *Invest Ophthalmol Vis Sci.* 2013 Jun 6;54(6):3912-21.
25. Yang J, Huang H, Yang H, He X, Jiang X, Shi Y, Alatangaole D, Shi L, Zhou N. Specific activation of the G protein-coupled receptor BNGR-A21 by the neuropeptide corazonin from the silkworm, *Bombyx mori*, dually couples to the G(q) and G(s) signaling cascades. *J Biol Chem.* 2013 Apr 26;288(17):11662-75.
26. Zheng C, Guo Q, Wu Z, Sun L, Zhang Z, Li C, Zhang X. Amphiphilic glycopolymer nanoparticles as vehicles for nasal delivery of peptides and proteins. *Eur J Pharm Sci.* 2013 Jul 16;49(4):474-82.
27. Sun X, Liu X, Zhang Y, Kuang X, Lv B, Ge J. A simple and effective pressure culture system modified from a transwell cell culture system. *Biol Res.* 2013;46(1):47-52.
28. Cai J, Han Y, Ren H, Chen C, He D, Zhou L, Eisner GM, Asico LD, Jose PA, Zeng C. Extracellular vesicle-mediated transfer of donor genomic DNA to recipient cells is a novel mechanism for genetic influence between cells. *J Mol Cell Biol.* 2013 Aug;5(4):227-38.
29. Li W, Zhou M, Ren H, Hu HM, Lu L, Cao M, Wang LX. Tumor-derived autophagosomes (DRibbles) induce B cell activation in a TLR2-MyD88 dependent manner. *PLoS One.* 2013;8(1):e53564.
30. Liu L, Zhao L, Yang J, Wan X, Hu N, Yeh L, Joo SW, Qian S. Low - Voltage Pulsed Electric Field Sterilization on a Microfluidic Chip. *Electroanalysis.* 2013 May;25(5):1301-9.
31. Liu H, Peng H, Wu Y, Zhang C, Cai Y, Xu G, Li Q, Chen X, Ji J, Zhang Y, OuYang HW. The promotion of bone regeneration by nanofibrous hydroxyapatite/chitosan scaffolds by effects on integrin-BMP/Smad signaling pathway in BMSCs. *Biomaterials.* 2013 Jun;34(18):4404-17.
32. Li H, Feng Z, Tsang TC, Tang T, Jia X, He X, Pennington ME, Badowski MS, Liu AK, Chen D, Harris DT, Martinez J, Meade-Tollin LC. Fusion of HepG2 cells with mesenchymal stem cells increases cancer associated and malignant properties: an in vivo metastasis model. *Oncol Rep.* 2014 Aug;32(2):539-47.
33. Qian J, Wu C, Chen X, Li X, Ying G, Jin L, Ma Q, Li G, Shi Y, Zhang G, Zhou N. Differential requirements of arrestin-3 and clathrin for ligand-dependent and -independent internalization of human G protein-coupled receptor 40. *Cell Signal.* 2014 Nov;26(11):2412-23.
34. Deng X, Yang H, He X, Liao Y, Zheng C, Zhou Q, Zhu C, Zhang G, Gao J, Zhou N. Activation of Bombyx neuropeptide G protein-coupled receptor A4 via a G α i-dependent signaling pathway by direct interaction with neuropeptide F from silkworm, *Bombyx mori*. *Insect Biochem Mol Biol.* 2014 Feb;45:77-88.
35. Mi XG, Song ZB, Wu P, Zhang YW, Sun LG, Bao YL, Zhang Y, Zheng LH, Sun Y, Yu CL, Wu Y, Wang GN, Li YX. Alantolactone induces cell apoptosis partially through down-regulation of testes-specific protease 50 expression. *Toxicol Lett.* 2014 Jan 30;224(3):349-55.
36. Yang X, Wu L, Duan X, Cui L, Luo J, Li G. Adenovirus carrying gene encoding *Haliothis discus discus* sialic acid binding lectin induces cancer cell apoptosis. *Mar Drugs.* 2014 Jun 30;12(7):3994-4004.
37. Song K, Han Y, Zhang L, Liu G, Yang P, Cheng X, Bu L, Sheng H, Qu S. ATP Synthase β -Chain Overexpression in SR-BI Knockout Mice Increases HDL Uptake and Reduces Plasma HDL Level. *Int J Endocrinol.* 2014;2014:356432.
38. Xu N, Wang J, Zhang ZF, Pang DW, Wang HZ, Zhang ZL. Anisotropic cell-to-cell spread of vaccinia virus on microgrooved substrate. *Biomaterials.* 2014 Jun;35(19):5049-55.
39. Wang M, Wang L, Guo Y, Sun R, Yue F, Yi Q, Song L. The broad pattern recognition spectrum of the Toll-like receptor in mollusk *Zhikong scallop Chlamys farreri*. *Dev Comp Immunol.* 2015 Oct;52(2):192-201.
40. Zhao X, Liu L, Hegazy AM, Wang H, Li J, Zheng F, Zhou Y, Wang W, Li J, Liu X, Lin L. Mannose receptor mediated phagocytosis of bacteria in macrophages of blunt snout bream (*Megalobrama amblycephala*) in a Ca(2+)-dependent manner. *Fish Shellfish Immunol.* 2015 Apr;43(2):357-63.
41. Zhou Y, Zhou Z, Zhang W, Hu X, Wei H, Peng J, Jiang S. SIRT1 inhibits adipogenesis and promotes myogenic differentiation in C3H10T1/2 pluripotent cells by regulating Wnt signaling. *Cell Biosci.* 2015 Nov 12;5:61.
42. Cao H, Zhang Z, Zhao S, He X, Yu H, Yin Q, Zhang Z, Gu W, Chen L, Li Y. Hydrophobic interaction mediating self-assembled nanoparticles of succinobucol suppress lung metastasis of breast cancer by inhibition of VCAM-1 expression. *J Control Release.* 2015 May 10;205:162-71.

43. Cai WB, Yang HL, Zhang J, Yin JK, Yang YL, Yuan LJ, Zhang L, Duan YY. The Optimized Fabrication of Nanobubbles as Ultrasound Contrast Agents for Tumor Imaging. *Sci Rep.* 2015 Sep 3;5:13725.
44. Zhang C, Yuan H, Liu H, Chen X, Lu P, Zhu T, Yang L, Yin Z, Heng BC, Zhang Y, Ouyang H. Well-aligned chitosan-based ultrafine fibers committed teno-lineage differentiation of human induced pluripotent stem cells for Achilles tendon regeneration. *Biomaterials.* 2015;53:716-30.
45. Wang D, Zhao J, Gao Z, Chen N, Wen B, Lu W, Lei Z, Chen C, Liu Y, Feng J, Wang JH. Neurons in the barrel cortex turn into processing whisker and odor signals: a cellular mechanism for the storage and retrieval of associative signals. *Front Cell Neurosci.* 2015 Aug 21;9:320.
46. Gu MJ, Li KF, Zhang LX, Wang H, Liu LS, Zheng ZZ, Han NY, Yang ZJ, Fan TY. In vitro study of novel gadolinium-loaded liposomes guided by GBI-10 aptamer for promising tumor targeting and tumor diagnosis by magnetic resonance imaging. *Int J Nanomedicine.* 2015 Aug 17;10:5187-204.
47. Lei Z, Liu B, Wang JH. Reward memory relieves anxiety-related behavior through synaptic strengthening and protein kinase C in dentate gyrus. *Hippocampus.* 2016 Apr;26(4):502-16.
48. Mi XG, Song ZB, Sun LG, Bao YL, Yu CL, Wu Y, Li YX. Cardamonin inhibited cell viability and tumorigenesis partially through blockade of testes-specific protease 50-mediated nuclear factor-kappaB signaling pathway activation. *Int J Biochem Cell Biol.* 2016 Apr;73:63-71.
49. Wan XK, Yuan SL, Tao HX, Diao LP, Wang YC, Cao C, Liu CJ. The Upregulation of TRAF1 Induced by *Helicobacter pylori* Plays an Antiapoptotic Effect on the Infected Cells. *Helicobacter.* 2016 Dec;21(6):554-564.
50. Yan YJ, Wang TM, Liu W, Wu CW, Zhu AY, Chi CF, Lü ZM, Yang JW. Identification and Expression Profile of the Gonadotropin-Releasing Hormone Receptor in Common Chinese Cuttlefish, *Sepiella japonica*. *J Exp Zool A Ecol Genet Physiol.* 2016 Aug;325(7):453-66.
51. Feng W, Gu YF, Nie L, Guo DY, Xiang LX, Shao JZ. Characterization of SIGIRR/IL-1R8 Homolog from Zebrafish Provides New Insights into Its Inhibitory Role in Hepatic Inflammation. *J Immunol.* 2016 Jul 1;197(1):151-67.
52. Tang Y, Cui Y, Li Z, Jiao Z, Zhang Y, He Y, Chen G, Zhou Q, Wang W, Zhou X, Luo J, Zhang S. Radiation-induced miR-208a increases the proliferation and radioresistance by targeting p21 in human lung cancer cells. *J Exp Clin Cancer Res.* 2016 Jan 12;35:7.
53. Lü ZM, Liu W, Liu LQ, Wang TM, Shi HL, Ping HL, Chi CF, Yang JW, Wu CW. Cloning, Characterization, and Expression Profile of Estrogen Receptor in Common Chinese Cuttlefish, *Sepiella japonica*. *J Exp Zool A Ecol Genet Physiol.* 2016 Mar;325(3):181-93.
54. Xu XG, Hu JF, Ma JX, Nie L, Shao T, Xiang LX, Shao JZ. Essential Roles of TIM-1 and TIM-4 Homologs in Adaptive Humoral Immunity in a Zebrafish Model. *J Immunol.* 2016 Feb 15;196(4):1686-99.
55. Li J, Gao J, Han L, Zhang Y, Guan W, Zhou L, Yu Y, Han W. Development of a membrane-anchored ligand and receptor yeast two-hybrid system for ligand-receptor interaction identification. *Sci Rep.* 2016 Oct 20;6:35631.
56. Fan X, Guo Y, Wang L, Xiong X, Zhu L, Fang K. Diagnosis of prostate cancer using anti-PSMA aptamer A10-3.2-oriented lipid nanobubbles. *Int J Nanomedicine.* 2016 Aug 12;11:3939-50.
57. Yang Z, Xiong F, Wang Y, Gong W, Huang J, Chen C, Liu P, Huang H. TGR5 activation suppressed S1P/S1P2 signaling and resisted high glucose-induced fibrosis in glomerular mesangial cells. *Pharmacol Res.* 2016 Sep;111:226-36.
58. Xie Y, Zhong Y, Gao T, Zhang X, Li LI, Ruan H, Li D. Human lymphatic endothelial cells contribute to epithelial ovarian carcinoma metastasis by promoting lymphangiogenesis and tumor cell invasion. *Exp Ther Med.* 2016 May;11(5):1587-1594.
59. Yang SD, Zhu WJ, Zhu QL, Chen WL, Ren ZX, Li F, Yuan ZQ, Li JZ, Liu Y, Zhou XF, Liu C, Zhang XN. Binary-copolymer system based on low-density lipoprotein-coupled N-succinyl chitosan liposomes for co-delivery of MDR1 siRNA and paclitaxel, enhances antitumor effects via reducing drug. *J Biomed Mater Res B Appl Biomater.* 2016 Mar 23. doi: 10.1002/jbm.b.33636.
60. Xu H, Li Z, Yu Y, Sizzdahkhani S, Ho WS, Yin F, Wang L, Zhu G, Zhang M, Jiang L, Zhuang Z, Qin J. A dynamic in vivo-like organotypic blood-brain barrier model to probe metastatic brain tumors. *Sci Rep.* 2016 Nov 10;6:36670.
61. Li W, Zhang L, Zhu Y, Zhang J, Shen L, Huang S, Fang S. Improving ovarian cancer imaging with LHRH-NBs: an experimental study. *Arch Gynecol Obstet.* 2016 Sep;294(3):631-8.
62. Li G, Gao Y, Cui L, Wu L, Yang X, Chen J. *Anguilla japonica* lectin I delivery through adenovirus vector induces apoptotic cancer cell death through interaction with PRMT5. *J Gene Med.* 2016 Apr; 18(4-6): 65-74.
63. Zhang X, Wang X, Zhong W, Ren X, Sha X, Fang X. Matrix metalloproteinases-2/9-sensitive peptide-conjugated polymer micelles for site-specific release of drugs and enhancing tumor accumulation: preparation and in vitro and in vivo evaluation. *Int J Nanomedicine.* 2016 Apr 22;11:1643-61.
64. Wang T, Yang Z, Zhou N, Sun L, Lv Z, Wu C. Identification and functional characterization of 5-HT4 receptor in sea cucumber *Apostichopus japonicus* (Selenka). *Sci Rep.* 2017 Jan 6;7:40247.

Version 2017.12.12