

DNA Ladders

Omega Biotek offers a wide range of products for the accurate size and mass determination of DNA. Most of DNA Ladders includes one or two higher-intensity bands for easy identification and orientation. The ladders are supplied premixed with loading buffer and are stable at room temperature for a few weeks. There is no need to heat, mix, or dilute ladders prior to loading them on gel.

1kb DNA Ladder

- **CAT.No.:** M03-01
- **Packing:** 250ul
- **Concentration:** 0.064 ug/ul
- **Store:** at 4°C for periods up to 6 months. For longer periods store at -20°C.
- **Description:** 1kb DNA Ladder is designed for sizing and approximate quantification of high range linear double-stranded DNA fragments on agarose gel. The ladder is composed of eight purified individual DNA fragments (in base pairs): 10000, 8000, 6000, 5000, 4000, 3000, 2000, 1000.
- **Direction:**
 - ◇ This product can load directly, since it contains 1× loading buffer. It is free of fluorescence dye such as EtBr.
 - ◇ For a loading well less than 6mm in width, 5ul of this product is enough for clear bands. For a wider well, more product may be required.
 - ◇ Gel and buffer may effect the electrophoresis. For better result, please use high quality agarose and fresh running buffer. 0.8% agarose gel is recommended.
- **Electrophoresis Map:**

